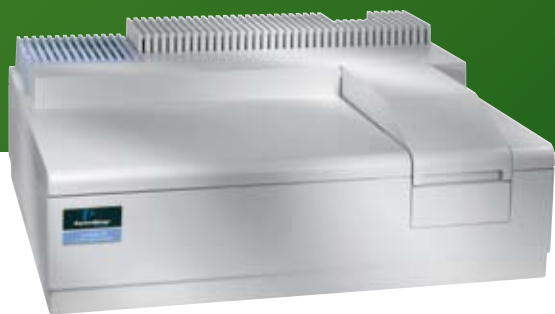
HUMAN HEALTH

ENVIRONMENTAL HEALTH

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8+1 Cell Changer installed in Lambda 25's sample compartment.

9+1 Cell Changer installed in Lambda 25's sample compartment.

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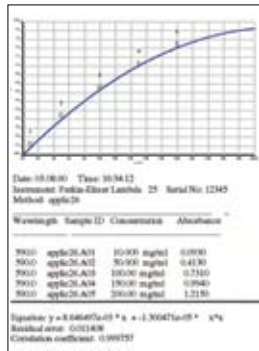
Table 1. Alternative cells for DNA analyses.

Cell Type	Part Number	Pathlength	Minimum Volume ¹	Nominal Volume
Semi-micro	80631064	10 mm	400 μL	1400 μL
Micro	80631077	10 mm	200 μL	400 μL
Ultra-micro	80631078	10 mm	50 μL	30 μL
Ultra-micro	80631061	10 mm	10 μL	5 μL
Ultra-micro	80631063	1 mm	10 μL	5 μL
Ultra-micro	80631060	5 mm	5 μL	2.5 μL
Ultra-micro	80631062	0.1 mm	2 μL	0.5 μL

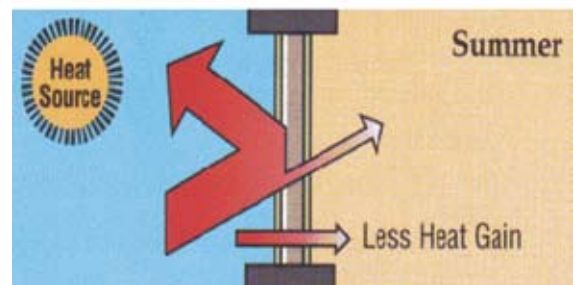
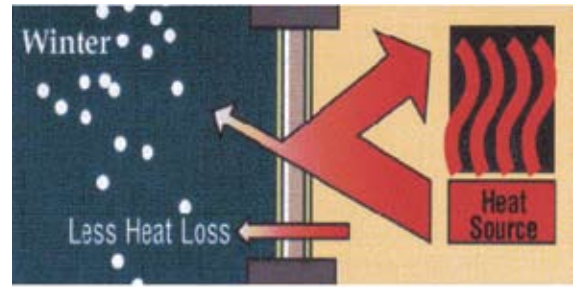
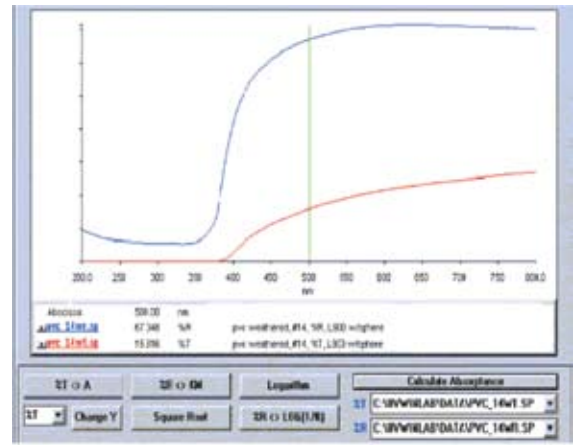
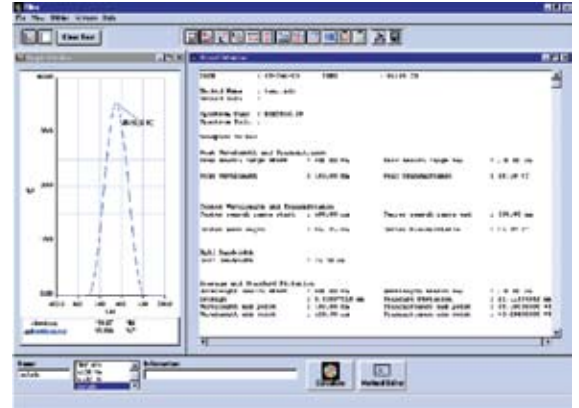
¹ The micro cells and ultra-micro cells must be "over-filled", so that the meniscus of the liquid is not in the sample beam. Therefore the minimum volume is larger than the "nominal volume". This problem does not apply to "normal" cells.

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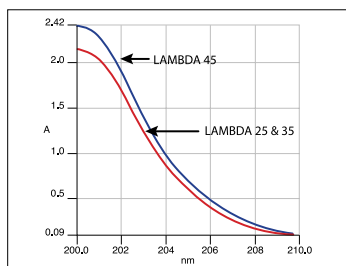
Figure 1. LAMBDA systems exceed the pharmacopeia stray light test specification of an absorbance of > 2.0A at 200 nm using a 1.2 w/v potassium chloride solution. The additional pre-monochromator in the LAMBDA 45 further reduces stray light compared to the LAMBDA 25 and 35.

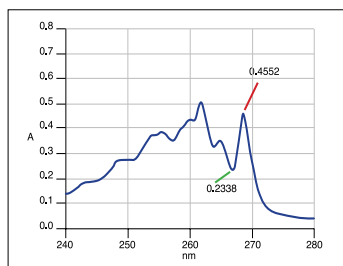
Figure 2. Pharmacopeia resolution test using 0.02% w/v toluene in hexane at 1 nm slit width. Ratio of peak and trough near 269 nm and 266 nm > 1.9 vs. the pass criterion of 1.5.

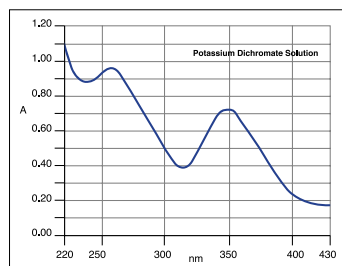
Figure 3. LAMBDA systems easily pass the pharmacopeia absorbance accuracy test at 235 nm, 257 nm, 313 nm and 350 nm using 60.06 mg/L potassium dichromate solution.

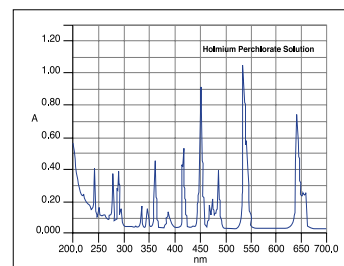
Figure 4. LAMBDA systems exceed pharmacopeia and ASTM requirements for wavelength accuracy (NIST 260-140) using a 4 mg/mL solution of holmium perchlorate.

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Copyright ©2011, PerkinElmer, Inc. 保留所有权利。PerkinElmer® 是 PerkinElmer, Inc. 的注册商标。