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made  
in  
Germany

**EMCLAB Instruments GmbH**

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“PC” means with PC software, Standard central beam height: 15 mm

M. Wave Professional PC-Control Software included in: EMC-16PC-V, EMC-16PC-UV

UV/VIS Analyst Software included in: EMC-31PC-UV, EMC-31PCS-UV  
EMC-61PC-UV, EMC-61PCS-UV

EMCLAB Spectrophotometers are tested with DAkkS\* certified UV/VIS Reference Materials  
NIST traceable:

- Photometric Accuracy 测光精度
- Wavelength Accuracy 波长准确度
- Spectral Resolution 光谱分辨率
- Stray Light 杂散光

\*Accreditation body for the Federal Republic of Germany

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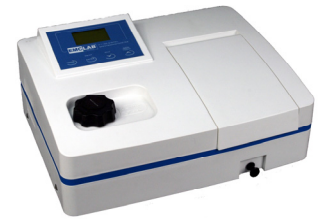
## UV / VIS Spectrophotometer EMC-16 series

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# UV / VIS Spectrophotometer EMC-16 series

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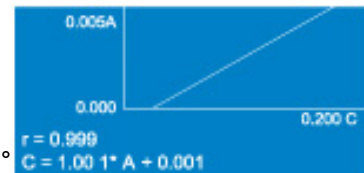
<input checked="" type="radio"/> Utility	
<input checked="" type="radio"/> D2 Lamp	On/Off
<input type="radio"/> W Lamp	On/Off
<input type="radio"/> Printer	

## UV / VIS Spectrophotometer EMC-3 series

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## UV / VIS Spectrophotometer EMC-6 series

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# NANO Spectrophotometer EMC-NANO-UV

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# UV / VIS Spectrophotometer EMC-3 & EMC-6 series

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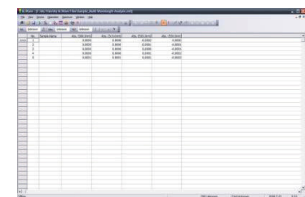
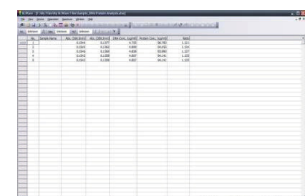
You can also enter other wavelengths and factors to calculate.

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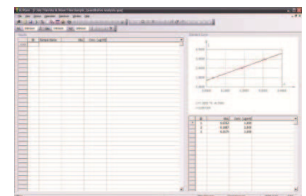
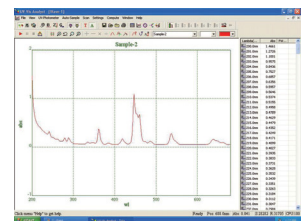
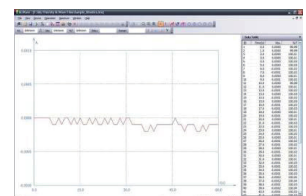
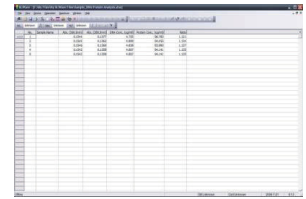
### DNA/Protein Test DNA/蛋白测试

Concentration and DNA purity are quickly and easily calculated. Absorbance ratios 260 nm/280 nm with optional subtracted absorbance at 320 nm

DNA concentration =  $62.9 \cdot A_{260} - 36.0 \cdot A_{280}$

Protein concentration =  $1552 \cdot A_{260} - 757.3 \cdot A_{280}$

Other wavelengths and factors may be entered.


Wavelength (nm)	Absorbance	Concentration
260	0.100	0.100
280	0.050	0.050
320	0.010	0.010

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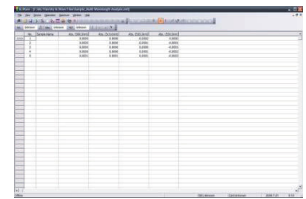
(auto cell changer is required to run multiple samples automatically)

Wavelength (nm)	Absorbance	Concentration
260	0.100	0.100
280	0.050	0.050
320	0.010	0.010

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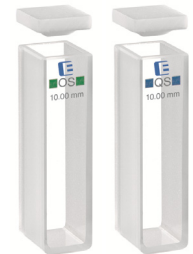比色皿	Standard	Premium
<b>Square</b>		
Glass 10 mm	EMC-S0001	EMC-100OS
Glass 50 mm	EMC-S0002	EMC-105OS
Quartz glass 10 mm	EMC-S0003	EMC-100QS
Quartz glass 50 mm	EMC-S0004	EMC-105QS
<b>Micro</b>		
Semi-micro cell glass 1000 $\mu$ l	EMC-S0005	EMC-108OS
Semi-micro cell quartz glass 1000 $\mu$ l	n/a	EMC-108QS
Semi-micro cell quartz glass 1000 $\mu$ l (black side walls and base)	n/a	EMC-108QSB
Micro cell quartz glass 500 $\mu$ l	n/a	EMC-108QSM
Micro cell quartz glass 500 $\mu$ l (black side walls and base)	n/a	EMC-108QSB
Ultra-micro cell quartz glass 100 $\mu$ l	EMC-S0008	EMC-108QSM15
<b>Flow-through</b>		
Flow-through cell glass 10 mm	EMC-S0009	EMC-175OS
Flow-through cell quartz glass 10 mm	EMC-S0010	EMC-175QS

## Spectrophotometer Accessories

### Cell Holder

4-cell holder standard for 10 mm square cuvette	EMC-00020
4-cell holder for up to 50 mm square cuvette	EMC-00021
Cell holder for micro cells, central beam height: 15 mm	EMC-00022
Single cell holder	EMC-00023
Single cell holder, water-jacketed	EMC-20024
4-cell holder, water-jacketed	EMC-00025

### Peltier Systems

Peltier System, Constant-Temperature-System, holder beam height: 15 mm	EMC-00026
Sipper System, holder beam height: 15 mm	EMC-00027
Peltier /Sipper System, Constant-Temperature System, holder beam height: 15 mm	EMC-00028

### Lamps

Halogen lamp 6V10W (only for EMC-11D-V)	EMC-00011
Halogen lamp 12V20W	EMC-00012
Halogen lamp 12V20W (only for EMC-6 series)	EMC-00013
Deuterium lamp for UV	EMC-00017

### Cleaner

Cleaning concentrate for cells/cuvettes and optical components, 1 L	EMC-S0020
---	-----------

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