LAC

Laboratory Furnaces
and Dryers
实验室电炉和干燥器

Art of heating

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S Dryers devices are designed for drying, vulcanization, preheating, curing and other low-temperature applications for treatment of various materials. The dryer is equipped with a chemically and mechanically-resistant stainless steel muffle to ensure long furnace life, and also boasts horizontal circulation of the internal atmosphere to provide uniform temperature distribution.

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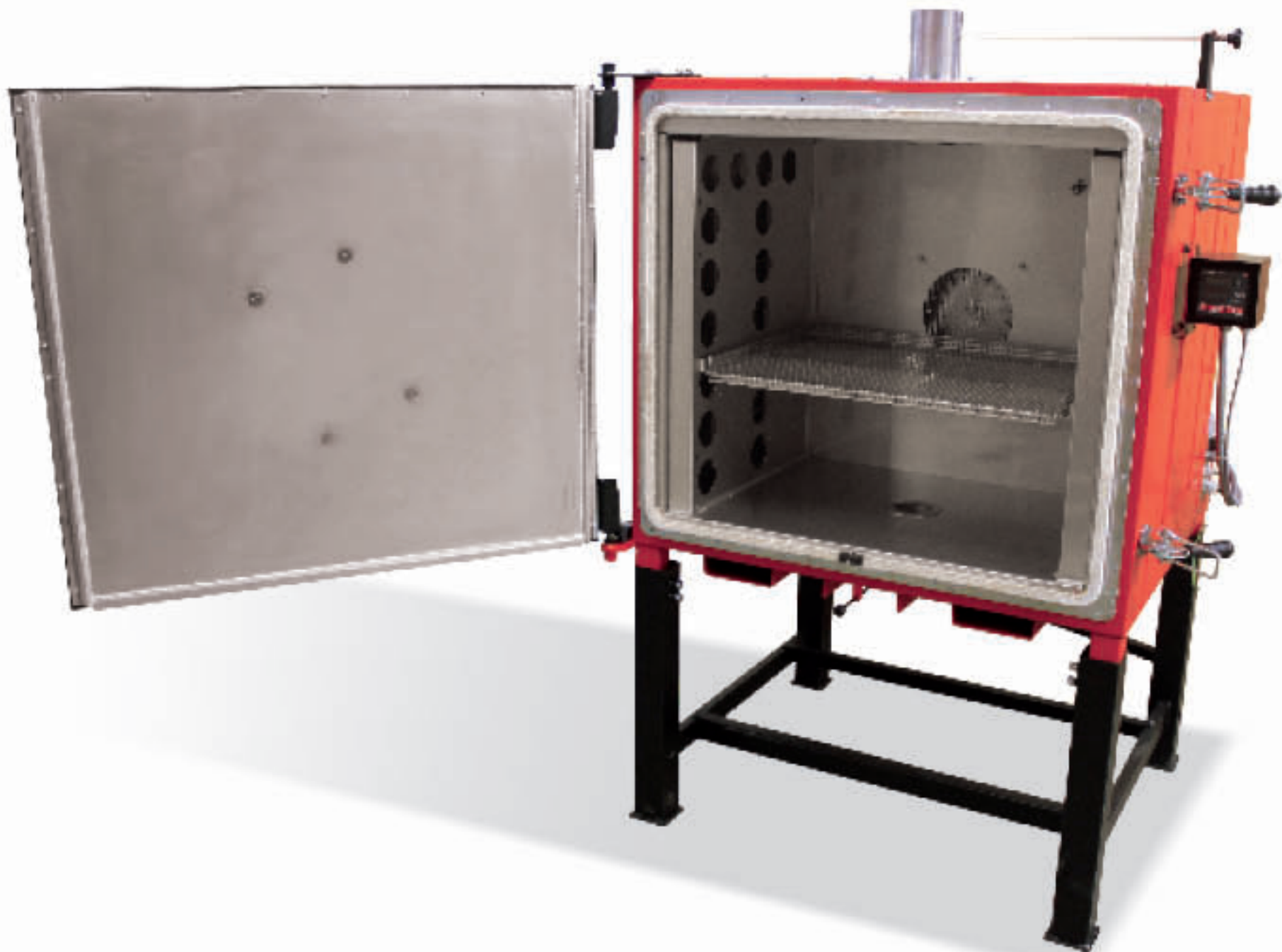
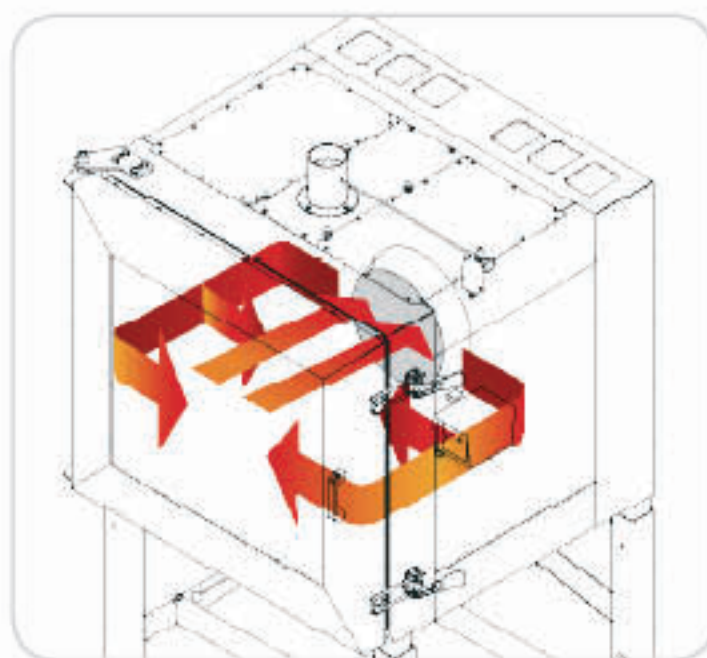
One of the greatest advantages of the PP tempering furnace is its horizontal air circulation of the inner atmosphere to ensure even heating of the batch. Uniform temperature distribution, precise regulation of temperature increase and decrease, and option of controlled cooling (for an additional fee) give the operator complete control of the process inside the furnace chamber. The PP furnace is robust and built to last for use in industry, but its compact space-saving size makes it also ideal for laboratory use. Treated with just a little care, your LAC PP furnace will serve you well for many years to come. This industrial furnace is tough enough for any lab.

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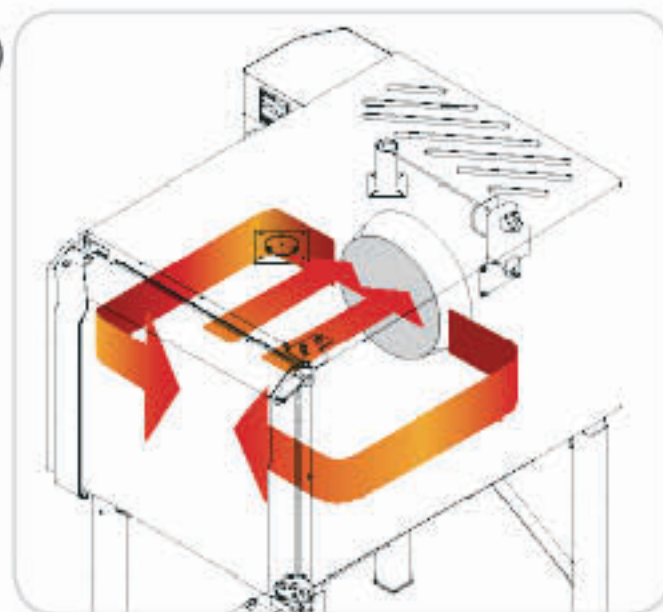
3D model of inner atmosphere circulation in PP furnaces

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Standard furnace equipped with:

- HtIndustry/Ht205 controller (30 programs)
- heating coils wound on ceramic tubes
- door sealing
- manually controlled ventilation flap
- manual door opening to the side
- limit unit
- inlet of protective atmosphere for one gas
- vacuum pressure gauge for vacuum control
- vacuum pump connection inlet
- thermistor for temperature control flange
- pressure relief valve with pressure 4-6 mbar
- ammeters to check heating element status
- type 'S' thermocouple
- silent contactless solid state relays

Precisely defined atmosphere protection to protect the surface of the charge from unwanted alteration and to ensure that the surface changes on the processed material are exactly as you require and nothing else. This furnace boasts low gas consumption; a great advantage when working with expensive gases like Argon, and of course the PKR can be set up for other standard gases such as nitrogen, forming gas, etc. If you would like to work at higher temperatures, an Inconel retort is available for an additional fee. The standard PKR refractory stainless retort can be used at temperatures up to 950 °C, the Inconel retort allows the furnace to be used at up to 150 °C more.

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GAS-TIGHT CHAMBER FURNACE WITH AIR CIRCULATION PKRC

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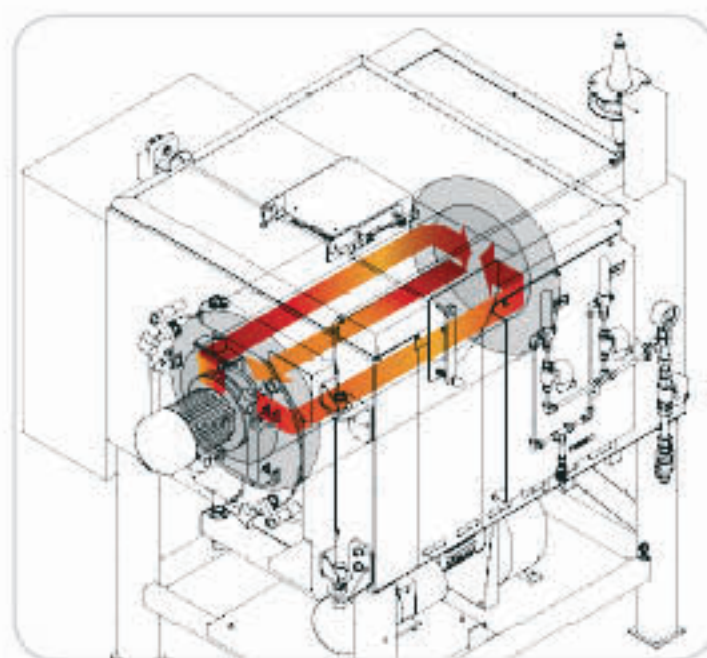
3D model of inner atmosphere circulation in PKRC furnaces

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The same applies to this furnace as to the PKR; it features precisely defined atmosphere protection to protect the surface of the charge from unwanted effects, low gas consumption, options for argon, nitrogen, forming gas, etc. In addition, the PKRC's circulation of the inner atmosphere gives it even better temperature distribution properties than the PKR. An Inconel retort is also available for the PKRC for an additional fee. A retort of this material allows the furnace to work at temperatures up to 150 °C higher than a standard stainless retort. The PKRC can also be adjusted to work as a nitriding furnace upon request.

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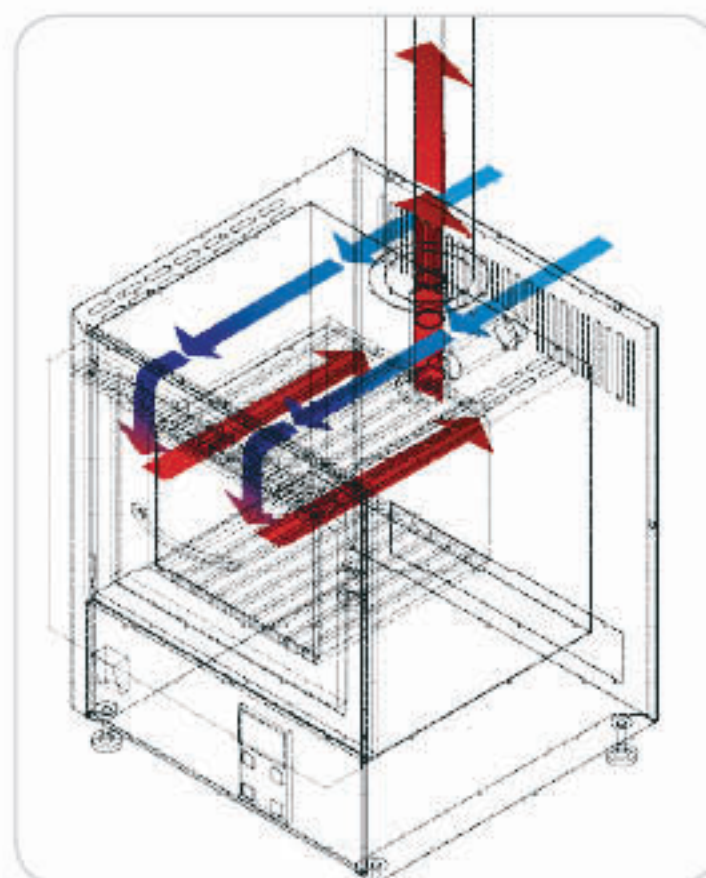
3D model of inner atmosphere circulation in LZ furnaces

Standard furnace equipped with:

- Ht40P controller (10 programs, 15 steps)
- heating elements in ceramic panels
- mineral fiber insulation panels
- type 'S' thermocouple
- contactless solid state relays for silent operation
- door limit switch for safe opening of the furnace

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- Ht205 controller, (30 programs, 15 steps)
- protective floor plate
- calibration of the measurement loop
- interface RS232 or EIA485 (including software)
- HTMonit set (software and interface)
- gas inlet for one or more inert gases

The LZ furnace ideal for the oxidation of samples and for testing to determine ash quantity. It is also suitable for any applications in which ample fresh air and also exhaust from the furnace atmosphere are required. Air is circulated throughout the internal furnace chamber and exchanged from six to twelve times per minute, depending upon the furnace temperature. The oxidation process is supported by the strong flow of oxygen. Air is preheated before it enters the LZ furnace chamber, thus allowing uniform distribution of heat even at high-temperature operation.

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Laboratories of all types appreciate the versatility of this furnace. Its heating elements are encased in ceramic panels, giving it extra protection from mechanical and chemical damage. L furnaces have a ventilation chimney flue for the extraction of fumes, but note that they are not suitable for pyrolysis. The design quality and materials used in the construction of the L furnace guarantee that it can handle even the most demanding tasks.

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Standard furnace equipped with:

- Ht40AL controller
(1 program: 2x ramp, 2x soak)
- muffle insulated with mineral fiber insulation mat
- heating coils wound around muffle, outside the chamber
- lid lined with insulation
- type 'S' thermocouple
- contactless solid state relays
- door limit switch for safe opening

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- HtIndustry / Ht205 controller (30 programs of 15 steps each)
- injector with exhaust fan and draft diverter for fume extraction
- inlet of protective atmosphere
- calibration of controller input measurement
- interface RS232 or EIA485 (including cable and software)
- HTMonit set (software and interface)
- gas inlet for one or more inert gases

This variation on the standard muffle furnace is for those who need to lower the charge into the furnace from the top. The heating coils are located outside of the furnace chamber where aggressive fumes cannot penetrate, and are protected by a tough ceramic muffle. The LMV furnace is also equipped with ventilation to protect from condensation during firing.

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LG 09/12

Standard furnace equipped with:

- Ht40P controller (10 programs, 15 steps)
- Digital scale with precision 0,01 g
- heating elements in ceramic panels
- ventilation chimney
- contactless solid state relays
- door limit switch for safe opening

Optional accessories (additional charge):

- Ht205 controller (15 steps, 30 programs)
- injector, exhaust fan and draft diverter
- inlet of protective atmosphere
- calibration of the measurement loop
- interface RS232 or EIA485
- HTMonit set (software and interface)

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The LG gravimetric furnace is a combination laboratory furnace, precision laboratory scale and software for the identification and evaluation of decreases in the charge weight. Measurements are output as a thermogravimetric curve documenting weight loss of samples as correlating to temperature and time. The LG furnace can test different biomass as well as certain types of waste such as RDF, plastics and paper etc.

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This hardening chamber furnace is right for you if you need to work at higher temperatures with higher volume batches. The PKE boasts speed and good heating accumulation into the brick lining, allowing you to open the furnace and check your charge more often. The furnace comes in desktop or freestanding designs, but don't let that fool you. The PKE is a tough furnace that brings industrial strength into your lab.

型号	最高温度	容量	外部尺寸 (wxdxh)	内部尺寸 (wxdxh)	输入	保护	重量*	电压	最大载重
	°C	l	mm	mm	kW	kg		V	kg
PKE 12/12	1280	12.2	700x650x850	250x200x250	3.0	95	16/1	230	20
PKE 18/12	1280	17.5	700x650x900	250x200x350	3.5	115	16/1	230	30
PKE 18/12R	1280	17.5	700x650x900	250x200x350	5.5	115	16/3	400	30
PKE 25/12	1280	25	700x1300**x1100	250x200x500	7.0	132	16/3	400	50
PKE 45/12	1280	44	800x1350**x1100	350x250x500	13.0	160	25/3	400	100
PKE 65/12	1280	65	800x1350**x1300	350x250x750	16.0	195	32/3	400	130
PKE 90/12	1280	87	800x1350**x1550	350x250x1000	18.0	225	32/3	400	150

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Standard furnace equipped with:

- Ht40AL controller (2x ramp, 2x soak)
- heating elements in insulation grooves
- mineral fiber insulation mats
- wiring located in a separate cabinet
- type 'S' thermocouple
- silent contactless solid state relays
- door limit switch for safe opening
- handle opens the furnace upwards
- desktop design

Optional accessories for additional charge:

- Ht205 controller (30 programs, 15 steps)
- injector, exhaust fan and draft diverter
- stand for vertical furnace placement
- inlet of protective atmosphere
- mineral fiber tube seal
- metal plugs for gas-tight design
- 3-zone heating system for 750mm tube
- calibration of controller input measurement
- interface RS232 or EIA485 (incl. software)

The constant temperature in the ceramic tube makes this furnace a researcher's best friend for trials and testing of various types of small samples. The charge can also be manipulated during testing without a major decline in temperature. The greatest advantage of the LT furnace is that the tube ends can be sealed with metal plugs. The plugs give the furnace gaslight properties so that you can work with the charge in protective atmosphere.

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Standard furnace equipped with:

- Ht205 controller (30 programs, 15 steps)
- heating coils in grooves
- lid insulated with mineral fiber mats
- door limit switch, contactless solid state relays

Optional accessories for additional charge:

- multi-channel controller for 10pt temp. scanning
- calibration of controller input measurement
- interface RS232 or EIA485 (including software)
- HTMonit set (software and interface)

Don't expect to save space or some kind of special design, but if you need to test different samples and scan the temperature of each all at the same time, then this furnace is definitely the right one for you. The LSP gradient furnace means you don't need five different furnaces, and you don't even have to keep changing your furnace settings. You just choose the temperature for the first zone, and the furnace will automatically set the proportional temperature in the other zones. Though you cannot regulate the temperature in the remaining zones, the temperatures are continuously scanned. If the temperature ratio in the individual zones is not as you would like, it can be adjusted upon request. The LSP gradient furnace is perfect for long-term sample testing as its firebrick lining has very good temperature accumulation properties.

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LH 15/13

Standard furnace equipped with:

- Ht40AL controller (2x ramp, 2x soak)
- manual door opening upwards
- heating elements built into grooves
- refractory brick lining
- ventilation chimney
- type 'S' thermocouple
- silent contactless solid state relay
- door safety limit switch

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The LH is a flexible furnace with a wide range of uses in the lab. This is the furnace you get when you need to test materials at higher working temperatures. The heating elements built into grooves in the furnace lining ensure very quick heat transmission. However, the LH is not suitable for treatment of aggressive materials as fumes can attack the unprotected heating coils or the firebrick lining.

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LABORATORY FURNACE WITH SILT RODS LHS

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LHS 08/15

Standard furnace equipped with:

- Ht205 controller (30 programs, 15 steps)
- silit rods on the furnace sides
- mineral fiber insulation panels
- ventilator forced cooling of the lining
- type ,B' thermocouple
- silent contactless solid state relays
- door safety limit switch
- manual door opening upwards
- desktop design

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Silit rods are an inexpensive way to reach higher working temperatures. They boast lower maintenance, a longer life, and are more resistant to temperature shock than the MoSi2 heating elements that are used in the VP furnace. The high-density silit rod construction of the LHS furnace allows it very rapid temperature increase. It can reach 1400 °C in 40 minutes. The LHS is also lighter and less bulky than the VP furnace.

型号	最高温度	容量	外部尺寸 (wxhxd)	内部尺寸 (wxhxd)	输入	重量	保护*	电压	最大载重
	°C	l	mm	mm	kW	kg		V	kg
LHS 03/14	1400	5	440x565x600	120x120x210	4	50	16/3	400	2
LHS 08/14	1400	8,4	490x615x680	170x170x290	8,0	60	16/3	400	5
LHS 15/14	1400	14,4	550x670x710	220x220x310	13	100	25/3	400	8
LHS 03/15	1500	5	440x565x600	120x120x210	4	50	16/3	400	2
LHS 08/15	1500	8,4	490x615x680	170x170x290	8,0	60	16/3	400	5
LHS 15/15	1500	14,4	550x670x710	220x220x310	13	100	25/3	400	8

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VP 10/16

VP furnaces are specially designed for high-temperature applications, but should be avoided when working at temperatures lower than 1200 °C. Use at such low temperatures will cause the VP to lose its precise temperature uniformity and can even lead to irreparable damage of the heating elements.

型号	最高温度	容量	外部尺寸 (wxdxd)	内部尺寸 (wxdxd)	最小输入	最大输入	重量	保护 **	电压	最大载重
	°C	l	mm	mm	kW	kW	kg		V	kg
VP 02/16	1600	2	660x680x740	130x150x135	3	3	90	16/1	230	1
VP 04/16	1600	4	660x680x740	130x160x180	3	3.5	80	16/1	230	4
VP 10/16	1600	10	850x1595*x775	200x200x250	6	9	290	25/3	400	6
VP 20/16	1600	20	850x1595*x775	250x250x310	8	12	315	32/3	400	10
VP 70/16	1600	70	1100x1750*x1150	400x300x600	14	18	350	63/3	400	20
VP 02/17	1700	2	660x680x740	130x150x135	3	3	90	16/1	230	1
VP 04/17	1700	4	660x680x740	130x160x180	3	3.5	80	16/1	230	4
VP 10/17	1700	10	850x1595*x775	200x200x250	6	9	290	25/3	400	6
VP 20/17	1700	20	850x1595*x775	250x250x310	8	12	315	32/3	400	10
VP 70/17	1700	70	1100x1750*x1150	400x300x600	14	18	350	63/3	400	20
VP 02/18	1800	2	660x680x740	130x150x135	3	3	90	16/1	230	1
VP 04/18	1800	4	660x680x740	130x160x180	3	3.5	100	16/1	230	4
VP 10/18	1800	10	850x1595*x775	200x200x250	6	9	290	25/3	400	6
VP 20/18	1800	20	850x1595*x775	250x250x310	8	12	315	32/3	400	10

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The CHTZ is a specially-designed work station for the heat treatment and chemical heat treatment of metallic materials, and it is not just for the laboratory. It is also particularly suited for applications where special emphasis is placed upon the quality of the surface treatment. The charge being treated remains in protective atmosphere throughout the entire heat treatment process, even while being transported to the hardening bath. This prevents the formation of oxide scaling on the surface of the charge. The CHTZ is suitable for hardening, carburizing, carbonitriding, nitriding and tempering.

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HARDENING WORKPLACE - SKM LABORATORY TABLE

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Our engineers have designed this modification of the LH 30/12 furnace especially for laboratory calcination of titanium dioxide (TiO₂). The furnace features a rotating quartz glass retort into which the charge is placed and then loaded into the furnace. The quartz retort, retort loader, and table are included with the furnace.

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This special version of the L laboratory furnace tests the thermal conductivity of samples at a Czech university. This is a bell furnace that features table-top bottom loading and has a 900mm high loading sill. The furnace chamber can be manually adjusted to tilt up to 300 mm. All the furnace walls are heated-up.

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This industrial research center furnace tests innovative high-strength steel and manufacturing technologies for the production of rail vehicles. It is primarily tasked with the heating of samples and temperature monitoring. Its components include a protective gas supply container with a thermocouple to enable batch processing in a protective atmosphere of argon, nitrogen or forming gas.

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PRODUCTION PLANT - INDUSTRIAL FURNACES AND DRYERS

LAC s.r.o.

Topolová 933,

667 01 Židlochovice

Czech Republic

PRODUCTION PLANT - REFRACTORY CASTABLE SHAPES

LAC, s. r. o.

Drnholecká 522

671 67 Hrušovany nad Jevišovkou

Czech Republic

BUREAU VERITAS
Certification

Certificate

Awarded to

LAC, s.r.o.

Head office: Štefánikova 116, 654 61 Rajhrad
Czech Republic

This is a multi-site certificate, additional site details are listed in the appendix to this certificate

Bureau Veritas Certification Holding SAS – UK Branch certifies
that the Management System of the above organisation has been audited
and found to be in accordance with the requirements of the management system
standard detailed below:

Standard

ISO 9001:2015

Scope of supply

DEVELOPMENT, MANUFACTURING, SALES AND SERVICE
OF FURNACES AND DRYERS.
DEVELOPMENT, MANUFACTURING, SALES AND SERVICE
OF OTHER GENERAL-PURPOSE MACHINERY.
MANUFACTURING AND SALES OF REFRACTORY PRODUCTS.

Original certification date: 28th NOVEMBER 2013

Recertification cycle start date: 28th NOVEMBER 2016

Subject to the continued satisfactory operation of the organisation's Management System,
this certificate expires on: 27th NOVEMBER 2019

Certificate number: CZ006822-2

Version 2, Revision date: 26th FEBRUARY 2016

Signed on behalf of BVZCH SAS UK Branch

Certification body address: 66 Piccadilly Street, London W1 9HQ, United Kingdom
Local office: BUREAU VERITAS CZECH REPUBLIC, spol. s r.o., Obvodská 1, 140 01 Praha 4, Czech Republic

0008

Further clarification regarding the scope of this certificate and the applicability of the management system requirements may be obtained
by consulting the registration. To check this certificate's validity please call: +420 210 088 216

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