

# Model KBF-S ECO 1020 | Constant climate chambers With thermoelectric cooling

Thanks to thermoelectric cooling technology with patented heat dissipation, the new KBF-S ECO climatic chamber is one of the most energy-efficient constant climate chambers on the market. The thermoelectric cooling technology means the KBF-S ECO series is also very quiet during operation.

#### **BENEFITS**

- Safe thanks to standard-compliant testing according to ICH Q1A guideline, even with full load
- Reliable due to failsafe operation without compromise
- · Smart as it is highly compatible for adaptation to diverse installation conditions and specific customer requirements
- · Economical as energy consumption is minimal thanks to optimized thermoelectric cooling



Model KBF-S ECO 1020



Model KBF-S ECO 1020

## **MAIN FEATURES**

- Temperature range (max. 24 °C under ambient temperature): o °C to +70
- Humidity range: 10% to 80% RH
- APT.line™ preheating chamber technology
- Independent water supply via tank
- LCD to display temperature and humidity along with additional information and alarms
- Internal data logger, measured values can be read out in open format via USB
- Unit self-test for comprehensive status analysis
- Inner chamber made completely of stainless steel
- 2 stainless steel racks, including shelf carrier
- Access port with silicone plug Ø 30 mm, left

#### ORDERING INFORMATION

Interior volume [L]	Voltage	Option model	Version	ArtNo.
1020	200230 V 1~ ph 50/60 Hz	Standard	KBFSECO1020-230V	9020-0419
	100120 V 1~ ph 50/60 Hz	Standard	KBFSECO1020UL-120V	9020-0422




## **TECHNICAL DATA**

Description	KBFSECO1020-230V <sup>1</sup>	KBFSECO1020UL-120V1
Article Number	9020-0419	9020-0422
Performance Data Temperature		
Temperature range (max. 24 °C unter Umgebungstemperatur) [°C]	070	070
Temperature variation at 40 °C [± K]	0.5	0.5
Temperature fluctuation at 40 °C [± K]	0.1	0.1
Max. heat compensation at 40 °C [W]	400	400
Performance Data Climate		
Temperature range (max. 24 °C unter Umgebungstemperatur) [°C]	570	570
Temperature variation at 25 °C and 60 % RH [± K]	0.1	0.1
Temperature variation at 40 °C and 75 % RH [± K]	0.5	0.5
Temperature fluctuation at 25 °C and 60 % RH [± K]	0.1	0.1
Temperature fluctuation at 40 °C and 75 % RH [± K]	0.1	0.1
Humidity range [% RH]	1080	1080
Humidity fluctuation at 25 °C and 60 % RH	o,4 % RH	0,4 % RH
Humidity fluctuation at 40 °C and 75 % RH	o,4 % RH	0,4 % RH
Recovery time after 30 seconds door open at 25 °C and 60 % RH [min]	5	5
Recovery time after 30 seconds door open at 40 °C and 75 % RH [min]	15	15
Electrical data		
Rated Voltage [V]	200230	100120
Power frequency [Hz]	50/60	50/60
Nominal power [kW]	1.2	1.2
Unit fuse [A]	16	16
Phase (Nominal voltage)	1~	1~
Measures		
Interior volume [L]	1020	1020
Net weight of the unit (empty) [kg]	322	322
Permitted load [kg]	150	150
Load per rack [kg]	45	45
Wall clearance back [mm]	100	100
Wall clearance sidewise [mm]	300	300
Internal Dimensions		
Width [mm]	976	976
Height [mm]	1250	1250
Depth [mm]	836	836
Doors		
Unit doors	2	2
Housing dimensions not incl. fittings and connections		
Width net [mm]	1254	1254
Height net [mm]	1926	1926
Depth net [mm]	1146	1146
Environment-specific data		
Energy consumption at 40 °C and 75 % RH [Wh/h]	135	135
Sound-pressure level [dB(A)]	49	49
Fixtures		
Number of shelves (std./max.)	2/15	2/15


<sup>1</sup> All technical data is specified for unloaded units with standard equipment at an ambient temperature of +22 °C ±3 °C and a power supply voltage fluctuation of ±10 %. The temperature data is determined to BINDER factory standard following DIN 12880, observing the recommended wall clearances of 10 % of the height, width, and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



# **DIMENSIONS** incl. fittings and connections [mm]

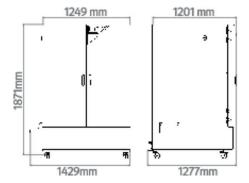

# **CHARTS**

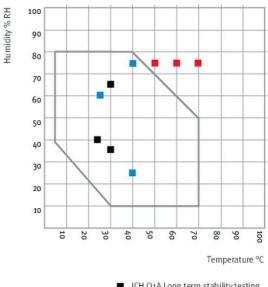

■ ICH Q1A Long term stability testing
■ ICH Q1A Accelerated stability testing

Climate chart

## **OPTIONS**

Designation	Description	*	ArtNo.
	left		
	30 mm	01	8012-2072
	50 mm	01	8012-2079
	100 mm	01	8012-2077
	right		
Access port with silicone	30 mm	01	8012-2071
lug	50 mm	01	8012-2078
	100 mm	01	8012-2076
	top		
	30 mm	01	8012-2073
	50 mm	01	8012-2074
	100 mm	01	8012-2075
Calibration certificate, expanded	for temperature and humidity; for extending the measurement in center of chamber to include another test value	-	8012-2066
Calibration certificate, temperature	temperature measurement incl. certificate and 27 measuring points at specified temperature	-	8012-2069
	temperature measurement incl. certificate, 15- 18 measuring points at specified temperature	-	8012-2068
	temperature measurement incl. certificate, 9 measuring points at specified temperature	-	8012-2067

<sup>\*</sup> Notes > See last page



Designation	Description	*	ArtNo.
Calibration certificate,	Measurement in center of chamber at 25 °C / 60% RH or at specified test values	-	8012-2065
temperature and humidity	temperature (according to DIN1288o) and humidity measurement incl. certificate, 27 temperature measuring points and 1 humidity measuring point, at 25 $^{\circ}$ C / 60 $^{\circ}$ RH or at specified values	_	8012-2070
Door lock	lockable door handle	-	8012-2087

<sup>\*</sup> Notes > See last page

# **ACCESSORIES**

Designation	Description	*	ArtNo.
APT-COM™ 4 BASIC-	for simple logging and documentation requirements with up to 5 networked units.		
dition	version 4, BASIC edition	-	9053-0039
APT-COM™ 4 GLP-	for working under GLP-compliant conditions. Measured values are documented in a tamper-proof way in line with the requirements of FDA Regulation 21 CFR 11.		
Edition	version 4, GLP edition	-	9053-0042
PT-COM™ <b>4</b>	convenient unit and user management built on the BASIC edition. Suitable for networking up to 100 units.		
ROFESSIONAL-Edition	version 4, PROFESSIONAL edition	_	9053-0040
SINDER PURE AQUA SERVICE	System for preparation or complete desalination of tap water, complete set containing PURE AQUA 300 single-use cartridge, measuring device, and all necessary connecting parts	_	8012-0759
SINDER PURE AQUA SERVICE, accessories	Single-use, replacement cartridge for BINDER PURE AQUA System	-	6011-0165
Convenient water supply	Flexible positioning of the water container with magnetic holder, hose adapter for water container with practical quick coupling, and water stop function		
	External, with magnetic holder	-	8012-1847
	T 220: For continuous temperature logging from -90 °C to 220 °C. The kit includes 1 data logger, Pt 100 sensor with 2 m extension cable and 1 magnetic fixture for mounting to the BINDER unit	-	8012-0715
Data Logger Kit	TH 70/70: For continuous temperature and humidity logging from -40 °C to 70 °C / 0% to 100% RH and additional logging of ambient conditions. Kit includes 1 data logger, 2 attachable combined humidity/temperature sensors, 2 m extension cable and 1 fixture for mounting to the BINDER unit	-	8012-0717
	TH 70: For continuous temperature and humidity logging from -40 °C to 70 °C; 0% to 100% RH. Kit includes 1 data logger, Pt 100 sensor with 2 m extension cable, and 1 magnetic fixture for mounting to the BINDER unit	_	8012-0716
ata Logger Software	LOG ANALYZE software kit, configuration and evaluation software for all BINDER Data Logger Kits (incl. USB data cable)	-	8012-082
	Basic set consisting of 2 pieces, attachment material, control unit for max. 4 light strips, 100-240 V, 50/60 Hz		
	Basic set 300, length 30 cm	-	8012-1107
	Basic set 500, length 50 cm	-	8012-1108
.ED light bars	Basic set 900, length 90 cm	-	8012-1109
LED light bais	Expansion set consisting of 2 pieces, attachment material: clips. For expanding the basic set of light bars		
	Expansion set 300, length 30 cm	-	8012-1716
	Expansion set 500, length 50 cm	-	8012-1717
	Expansion set 900, length 90 cm	-	8012-1718
H-neutral detergent	concentrated, for gentle remove of residual contaminants; 1 kg	-	1002-0016
	IQ/OQ documents – supporting documents for validation performed by customers, consisting of: IQ/OQ checklists incl. calibration guide and comprehensive unit documentation; parameters: temperature and humidity values		
	Digital in PDF format	-	7057-0002
Dualification documents	Hard copy inside folder	-	7007-000
qualification documents	IQ/OQ/PQ documents – supporting documents for validation performed by customers, according to customer requirements, PQ section added to qualification folder IQ/OQ; parameters: temperature and humidity values		
	Digital in PDF format	-	7057-000
	Hard copy inside folder	-	7007-000
Rack	stainless steel	-	6004-014
Rack accessories	fasteners (1 set of 4) for additional security of racks	-	8012-062
Rack, reinforced	stainless steel, with fasteners (1 set of 4)	-	8012-096
Shelf, perforated	Stainless steel	-	8009-079
WLAN kit	The kit contains a client bridge that establishes a wireless connection between BINDER units and APT-COM4, LIMS, or customer-specific software via an Ethernet interface. It provides an alternative solution in situations where units need to be positioned in locations without an on-site Ethernet connection. In secured networks, installation and configuration must be performed by the customer's IT service.	-	8500-0164

<sup>\*</sup> Notes > See last page



## **SERVICES**

Designation	Description	*	ArtNo.
Installation services			
Unit commissioning	Connect the unit to the customer-side connections (electricity, water, wastewater, gas), basic functions check, brief operating instructions. (excl.: unpacking, setup, controller instructions, programming, installation work)	05	DL10-0300
Unit instructions	Instruction regarding operating principle and basic functions of the unit, operation of the control electronics including programming	05	DL10-0700
Maintenance contracts			
Maintenance services			
BRONZE 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts	05	DL20-0710
GOLD 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts, testing of all key functions, replacement of wear parts, calibration of one temperature/humidity value, including certificate	05	DL20-0920
Maintenance	One-off maintenance service in accordance with maintenance schedule. Visual inspection of mechanical and electrical components, testing of all key functions. Calibration of a test temperature specified by the user in center of usable space without certificate	05	DL20-040
SILVER 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts, testing of all key functions, calibration of one test temperature specified by the user in the center of the usable space, without certificate	05	DL20-082
Calibration services			
	Expansion –Temperature and humidity calibration with 1 measuring point in center of chamber with 1 specified pair of values, including certificate	03, 04, 05	DL30-0302
Temperature and humidity calibration	Temperature and humidity calibration with 1 measuring point in center of chamber with 1 specified pair of values, including certificate	03, 04, 05	DL30-0301
Temperature and humidity measurement according to DIN12880	Temperature measurement in accordance with DIN 12880 with 27 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0427
Temperature and humidity measurement, 18-1 measuring points	Temperature measurement with 18 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0318
Temperature and humidity measurement, 27-1 measuring points	Temperature measurement with 27 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0327
Temperature and humidity measurement, 9-1 measuring points	Temperature measurement with 9 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0309
Validation services			
Execution of IQ/OQ	Execution of IQ/OQ in accordance with qualification folder	05	DL42-030
Execution of IQ/OQ/PQ	Execution of IQ/OQ/PQ in accordance with qualification folder	05	DL44-050
Warranty service			
1-year warranty extension	The warranty is extended by 1 year from the delivery date, wear parts are excluded	-	DL50-0030
+ NI=+== C== l==+ ====			

<sup>\*</sup> Notes > See last page



#### **NOTES**

- Condensation may occur in the area around the access port. Access ports may be placed in custom locations for an additional charge. UL mark is not granted when this option is used. 01
- 02
- Sensor calibration is performed in an accredited calibration laboratory. 03
- Calibration is performed according to the BINDER factory standard.
- Quoted prices do not include travel costs. Please refer to the chapter on BINDER Service for travel costs for your region. Quoted prices for services performed in Switzerland do not include a country-specific added fee (available on request).

## **BINDER GmbH**

Tuttlingen, Germany TEL +49 7462 2005 0 FAX +49 7462 2005 100 info@binder-world.com www.binder-world.com

BINDER Asia Pacific (Hong Kong) Ltd.

Kowloon, Hong Kong, P.R. China TEL +852 39070500 FAX +852 39070507 asia@binder-world.com www.binder-world.com

**BINDER Environmental Testing** Equipment (Shanghai) Co., Ltd. Shanghai, P.R. China TEL +86 21 685 808 25 FAX +86 21 685 808 29 china@binder-world.com www.binder-world.com

**Representative Office for CIS** 

Moscow, Russia TEL +7 495 988 15 16 FAX +7 495 988 15 17 russia@binder-world.com www.binder-world.com

**BINDER Inc.** 

Bohemia, NY, USA TEL +1 631 224 4340 FAX +1 631 224 4354 usa@binder-world.com www.binder-world.us