JUNIOR

Reaction Screening OSR

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UNCHAINED
LABS

Cover all screening and optimization bases 涵盖全部筛选和优化变量

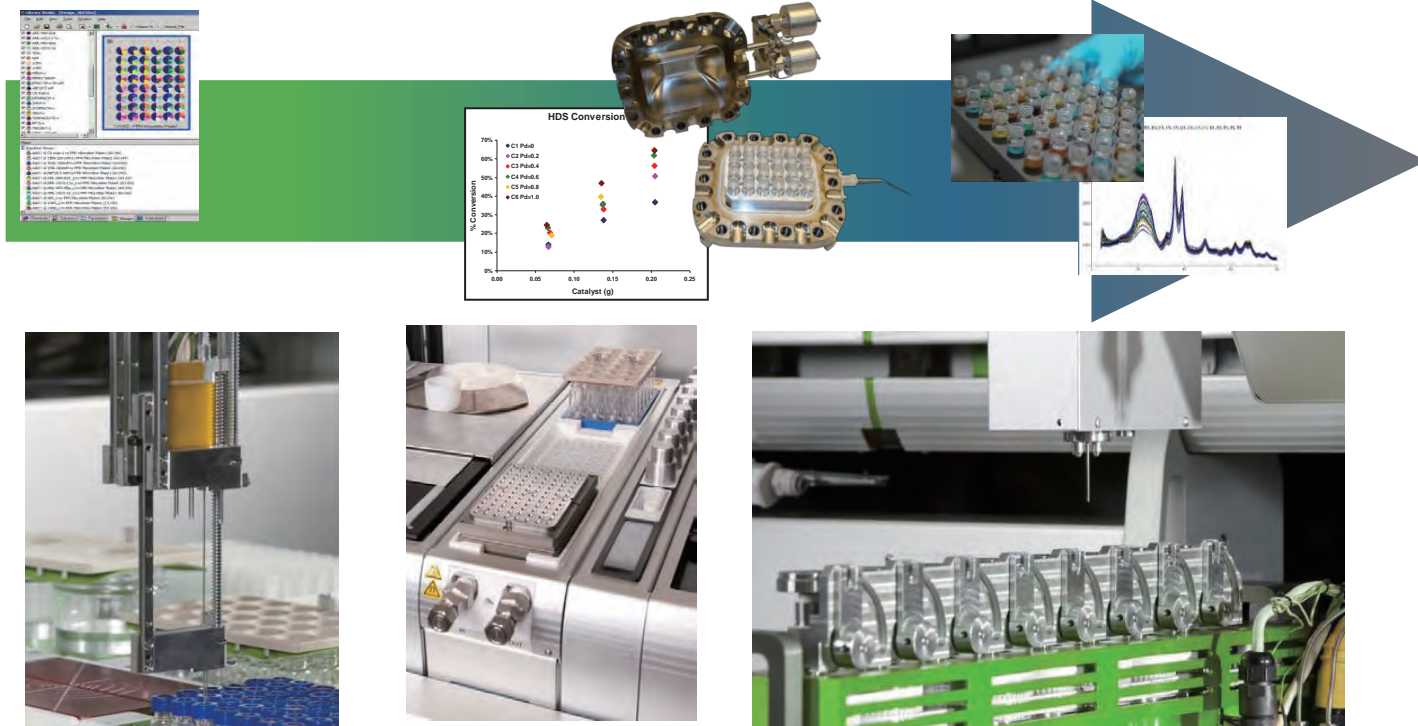
Junior screens all your substrates, catalysts, reactants, solvents and reaction conditions so you can tackle a huge range of organic transformations. Tight reaction control, stir plate temps, filtration and dilution mean your samples are ready for structural analysis whenever you are. Junior turns reaction screening and process optimization into totally routine activities. Build a software protocol for reaction screening and plow through hundreds of modifications per week, design real-time sampling to fine tune your chemistry pathways. Junior lets you dive deep so that you know exactly what tweaks to make next.

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Perform Multiple Applications 执行多种应用

Junior fully automated critical chemistry workflows to allow for you to walk up, set up your runs, and walk away worry-free.

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Reaction screening and optimization

Process optimization

Real-time reaction kinetics

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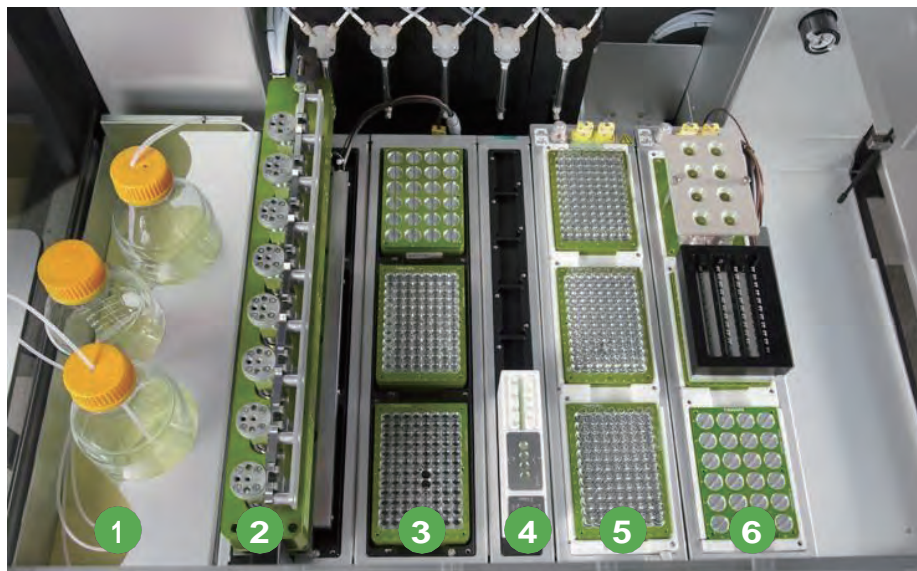
Dial in your Screening Process 了解您的化学反应

Junior lets you check out a ton of reaction screening variables like solvents, ligands, catalysts and reagents all at once. It'll plow through hundreds of modifications per week. Outfit one with an Optimization Sampling Reactor (OSR) to grab samples from your reactions as they're happening and know even more right now.

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

OSR Heated Tip High Pressure Sampling OSR高温高压取样针设计

Going the one variable at a time route doesn't cut it when you need to know how all your different reaction variables play together. Add an Optimization Sampling Reactor (OSR) to Junior and get real-time kinetics on all your reactions. OSR grabs time-point samples from up to 8 pressure and temp-controlled vessels at a time without interrupting a single reaction. Each vessel makes sure the heating, cooling and stirring for each of your reactions is just right.

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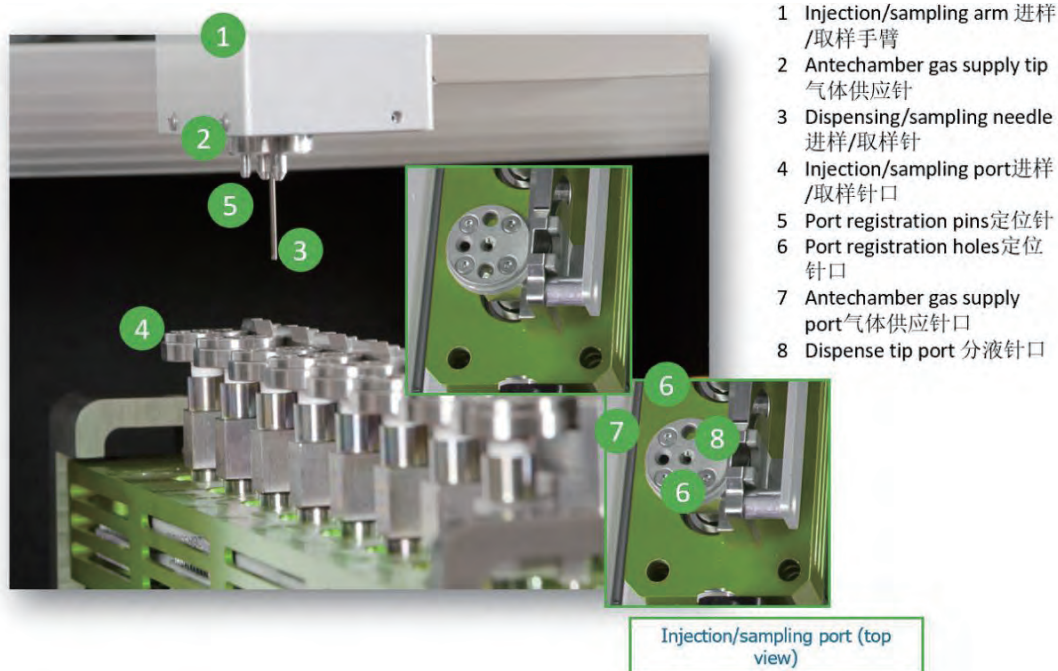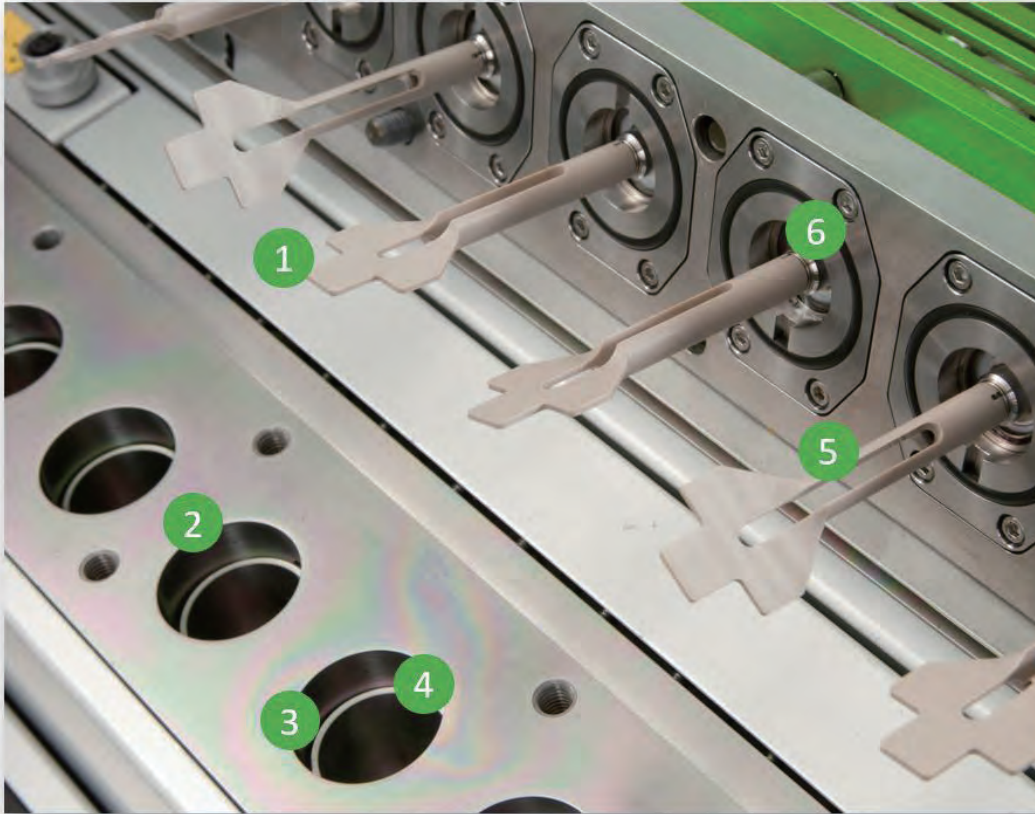
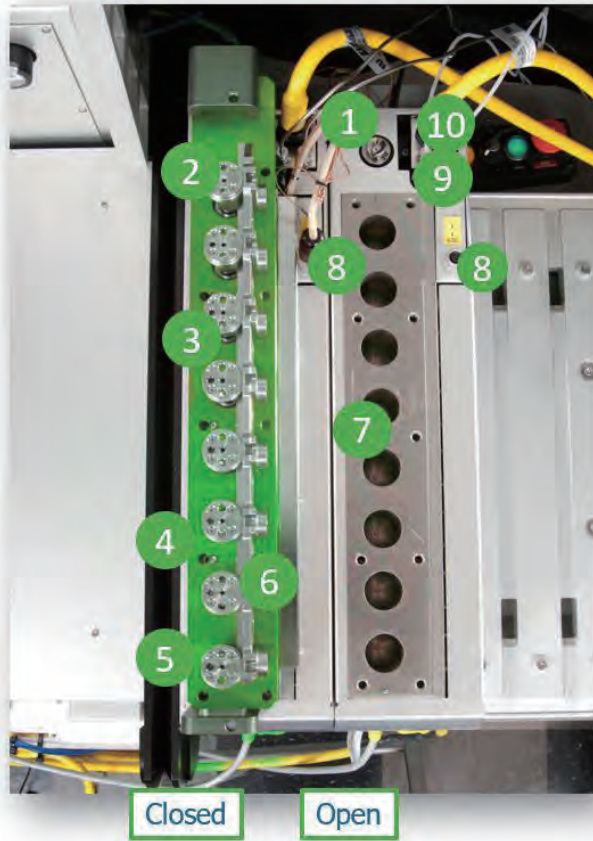
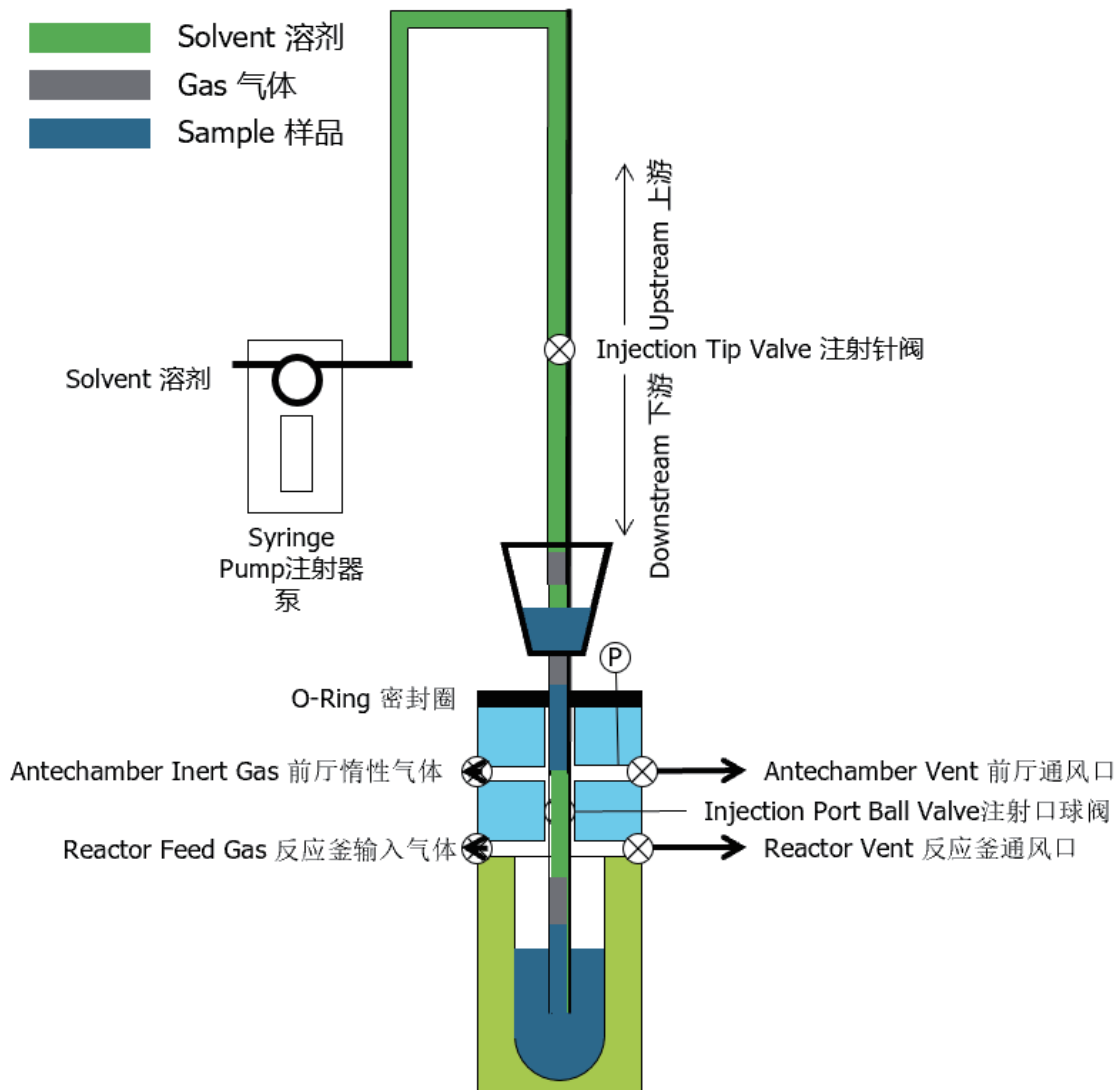
OSR In-situ Sampling Illustration OSR 高压取样图解


OSR In-situ Sampling Illustration OSR高压取样图解

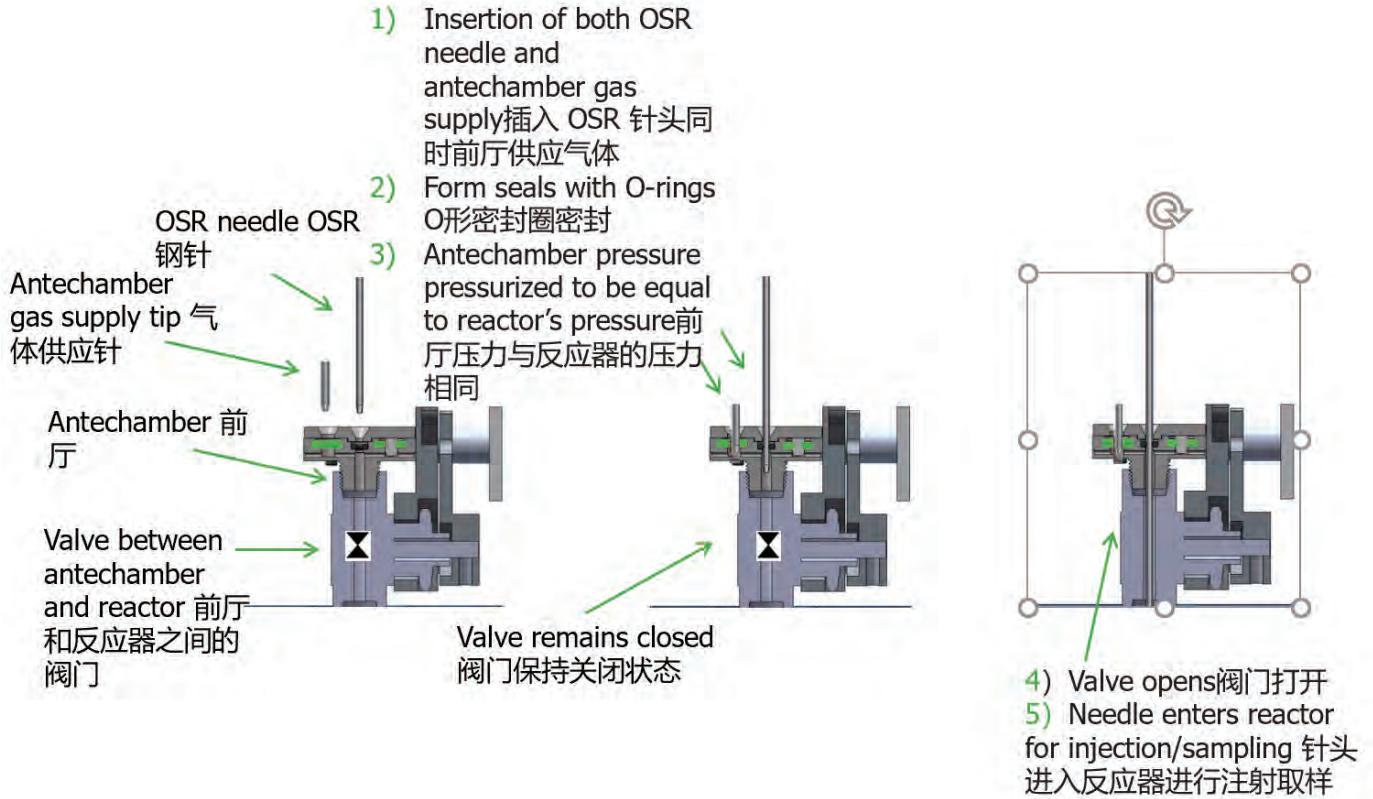Software break through bottlenecks 软件优势



LEA doesn't move the bottleneck, it totally unclogs it. You get intuitive experimental design that makes sense on the front end and full integration with analytical tools on the back end. LEA also links your conditions, steps and analytical data together, so your final report doesn't just have numbers, it's got all the information you need to make real decisions.

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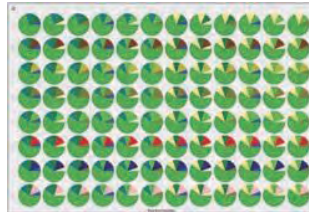Library Studio
Design complex, high-throughput experiments in an array-based format
设计复杂高通量实验

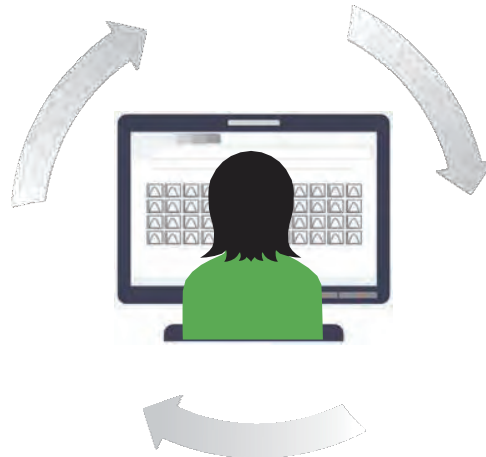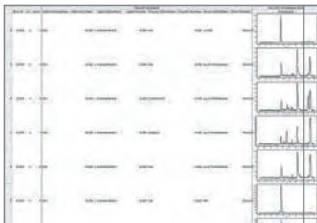PolyView
Review and report all info from experimental design, execution and analytics
查阅实验执行和结果的报告

Automation Studio
Execution of designed experiments and integrated analytics
执行实验和分析工具

Junior Reaction Screening OSR Specifications Junior 反应筛选仪参数

Heated tip liquid dispenser—extended tip

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Heated reservoir volume per tip: 1mL

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Reservoir temperature: Up to 120°C

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Temperature uniformity: $\pm 2^\circ\text{C}$ (across tips)

温度均匀性: $\pm 2^\circ\text{C}$

Tip pitch: 9mm

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Extendible tip: 1

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Syringe sizes: 50 μL –2.5 mL (standard supplied)

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Heated single-tip liquid dispenser

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Orbital: 60–3570rpm 转速

Maximum vortexing mass: 2268g(5lb/plate) 涡旋器能承受最大重量

Optimization sampling reactor and sampling arm 优化反应取样器和取样臂

Temperature rating: -20–200°C 温度范围

Reactors: 8 independent reactors 反应釜: 8 个独立反应釜

Possible delta between adjacent reactors: 50°C 相邻反应釜之间的可能温差: 50°C

Heating rate: Up to 5°C/min

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Cooling rate: Up to 2°C/min

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Reactor total volume: 40mL

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Reactor working volume: 5–25mL

Junior Reaction Screening OSR Specifications Junior 反应筛选仪参数

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Maximum overhead stirringspeed:750rpm

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Pressure rating:30–400psi

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Off-deck third-party instrument integration option

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Other systems available for virtual integration.

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Please contact Unchained Labs for a full list of systems.

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With integrated enclosure: 外壳尺寸

105cmWx90.4cmDx140cmH, ~150kg

With integrated table option: 工作台尺寸

167cm Wx 90,4 cm D x 200cm H, ~240 kg

Electrical:

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Junior:

120–220V±10%,50–60Hz,16A

Computer: 电脑

US:115V±10%,60Hz

EU:220V±10%,50Hz

Compressed dry air: 0.5MPa to 0.9MPa (70–130psi), 4L/min (8mm hose)

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