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Condenser ultrasound microphone Avisoft-Bioacoustics CM16/CPMA
Externally polarized condenser microphone for UltraSoundGate xx16x

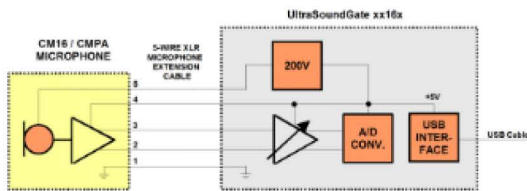
tripod


- 1, Frequency range : 2 kHz - 200 kHz
- 2 , Approximate input-referred self-noise Level (bandwidth: 30-50 kHz): 18 dB SPL
- 3, Approximate sensitivity: 500mV/Pa
- 4, Connector: male XLR-5 plug
- 5, Removeable protection grid
- 6, Dimensions : 36 x 36 x 60 mm
- 7, Weight : 65 g
- 8 , Supplied accessories: 5 m extension cable (other lengths of up to 300 m are available on request), high-quality 9, Sennheiser microphone clamp and mini
- 10, Optionally available USB-powered microphone heating

The condenser microphone Avisoft Bioacoustics CM16/CPMA provides a good compromise between sensitivity and flatness of the frequency response. It consists of an externally polarized metalized film diaphragm and a metal back plate. The capsule is replaceable in order to handle the relatively low durability of the thin diaphragm. This microphone model requires an additional 200V polarization voltage that is supplied by the UltraSoundGate base units. An optional USB-powered heating module can prevent condensation when the microphone is being used in the field.

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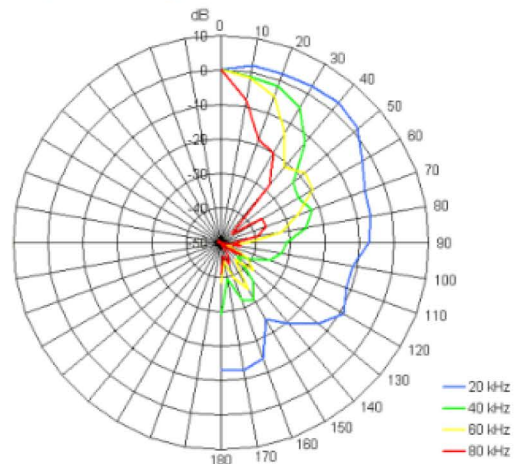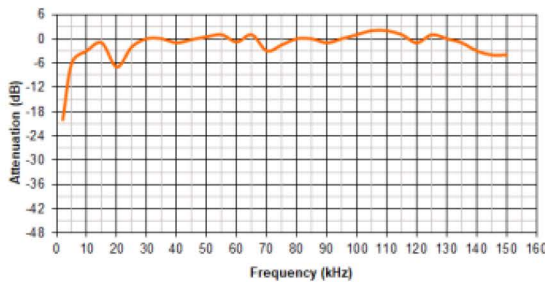
Avisoft Bioacoustics UltraSoundGate

Single-Channel Recording Systems

USG 116H : full-featured unit
sample rates of up to 1 MHz with external trigger inputs and spare microphone capsule (order # 51163 / ..4)

USG 116Hb : basic and cost-effective solution
sample rates of up to 300 kHz, without external trigger inputs and without spare microphone capsule (order # 51161 / ..2)

Special microphones for best results

The supplied sensitive externally polarized solid-dielectric condenser microphones exhibit a fairly flat frequency response from 15 to 200 kHz (order # 40011)

Alternative microphone models are available in order to suit special needs.

Functional principle of the single-channel system

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