

**EASY, SAFE AND REPRESENTATIVE
MEASUREMENTS WITH TIMEGATED® RAMAN**

SAMPLECUBE



EYE SAFETY

Raman techniques often use powerful laser sources which emphasizes the importance of proper eye safety measures. The SampleCube sample enclosure can be closed using a sliding door minimizing the risk of any high energy laser light causing retinal damage. The inside of the sample enclosure is colored black which reduces light reflections.

SAMPLING A LARGER SURFACE AREA

The SampleCube enables measurements from a larger sample surface for more representative measurement results. During measurements, the sample cup holder has an option for eccentric rotation and the measurement probe can be simultaneously moved in a scanning motion. This way a larger surface area is included in the sampling volume instead of a single point measurement. The user may choose to only sample a smaller surface area from the center of the sample or adjust the amplitude setting to include the whole sample surface area.

LIMITING LASER EXPOSURE

The scanning cycle is designed to cover the whole surface area effectively while limiting laser exposure of any single point. Because of this, many of the more sensitive samples can be measured, as sample degradation caused by laser exposure is minimized. The user may adjust the scanning speed to even further reduce the longest continuous time a single point is exposed to laser light.

EASY FOCUS DISTANCE ADJUSTMENT

The SampleCube also enables easy laser focusing with all compatible measurement probes. The probe can be moved up and down in a controlled manner using the adjustment screw on top. Once a suitable position has been found, the probe will stay in the chosen position during measurements.

SUITABLE FOR MANY SAMPLE TYPES

The main use case for the SampleCube are powder samples but it can also be used to measure other flat or flattenable samples. The SampleCube can also be used to measure liquid samples in vials or cuvettes as the measurement probe can be inserted horizontally in the sample enclosure. Using the SampleCube, all the aforementioned sample types can be measured in a controlled and eye safe manner.

ENHANCED EYE SAFETY

Easy Timegated® Raman measurements that cover a larger sample surface area while providing enhanced eye safety.



Closable chamber increases eye safety

MORE REPRESENTATIVE SAMPLING

The measured surface area can be adjusted for more representative sampling.



Easy laser focusing and probe adjustment

 **timegate**

Find us on:



@TgiRaman



Timegate Instruments Ltd

Visiting address:

Timegate Instruments Ltd

Tutkijantie 7

FI-90590 OULU, FINLAND

info@timegate.com

www.timegate.com



SAMPLECUBE

- A built-in display makes SampleCube adjustments easy
- Compatible with B&W TekBAC 100 series probes or any other probes with similar dimensions
- The scanning speed can be adjusted for improved sampling and decreased single point laser exposure
- **Time-resolved Raman analysis with SampleCube**

EASY, SAFE AND REPRESENTATIVE
MEASUREMENTS WITH TIMEGATED® RAMAN

 **timegate**



北京海菲尔格科技有限公司
联系人：孙经理
地址：北京市朝阳区南磨房路37号10层1008室
电话：010-53779530、13716489005
网址：www.hiferg.com
邮箱：nina.sun@hiferg.com

Visiting address:

Timegate Instruments Ltd
Tutkijantie 7
FI-90590 OULU, FINLAND
info@timegate.com
www.timegate.com