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## Controller screens are easy to navigate for fast and easy operation.

The AutoBlock can be operated in Manual Mode, allowing the analyst to perform a variety of functions independently.

The independent racking system enables the analyst to perform up to three digestion methods simultaneously.

Using the AutoBlock software, the analyst can easily create and save custom-designed methods.

The AutoBlock software will alert the analyst and perform a variety of maintenance and calibration procedures as required.

Accurately and efficiently add concentrated acids to 54 samples for accurate, hands-free digestions in a HEPA filtered environment.

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UNITED STATES ENVIRONMENTAL PROTECTION AGENCY  
WASHINGTON, D.C. 20460

MAR 31 1998

OFFICE OF  
WATER

John Stone  
Vice President  
Environmental Express  
490 Wando Park Blvd.  
Mt. Pleasant, SC 29464

Dear Mr. Stone:

This letter responds to your letter dated February 23, 1998 regarding the use of the Hot Block™ digestion system for EPA approved methods that use conventional hot plate digestion technique. EPA does not approve or endorse specific instrumentation. However, EPA does not restrict the use of instruments or equipment that employ technology which is different from the EPA approved methods as long as the user follows the procedures required in the EPA methods, as well as the manufacturer's operating instructions for proper use. Therefore, provided the above conditions are met, the Hot Block™ digestion system which is manufactured by your company is acceptable for use with EPA approved methods.

If I can be of additional assistance on this and other matters, please write or call me at (202) 260-1639 at your convenience.

Sincerely,

UNITED STATES ENVIRONMENTAL PROTECTION AGENCY  
WASHINGTON, D.C. 20460

MAR - 8 2001

OFFICE OF  
WATER

John Stone  
Vice President  
Environmental Express, Ltd.  
490 Wando Park Blvd.  
Mt. Pleasant, SC 29464

Dear Mr. Stone:

This letter responds to your letter regarding the use of the Environmental Express HotBlock™ digestion system with EPA methods currently approved or soon to be proposed for the determination of mercury. Specifically your letter referenced EPA Methods 245.1, 245.2, 245.7 (to be proposed), 7470, 7471, and 1631. As stated in an earlier letter, EPA does not approve or endorse specific instrumentation. However, provided that all procedures in the approved methods are followed, and all QA/QC criteria are met, the HotBlock™ is acceptable for use as a heating source in methods approved or proposed by EPA's Office of Water for use in National Pollutant Discharge Elimination System (NPDES) or National Primary Drinking Water Regulations (NPDWR) compliance monitoring. These methods include but are not limited to EPA Methods 245.1, 245.2, 245.7 and 1631.

For written clarification regarding the use of the Environmental Express HotBlock™ with EPA Methods 7470 and 7471, or other methods contained EPA's Office of Solid Waste methods manual, known as Test Methods for Evaluating Solid Waste: Physical/Chemical Methods (SW-846), you will need to contact EPA's Office of Solid Waste. Written inquiries regarding SW-846 methods may be submitted to the following address: OSW Methods Team, U.S. Environmental Protection Agency Office of Solid Waste (5307W), Ariel Rios Building, 1200 Pennsylvania Ave. NW, Washington, DC 20460, Fax: 703-308-0511.

Alternatively, you may submit an inquiry through the Methods Information Communication Exchange (MICE) Service. The service provides timely answers to method-related questions and takes comments on technical issues regarding the EPA Office of Solid Waste methods. The MICE Service can be reached by telephone (703-676-4690), by fax (703-903-1373), or by e-mail: [mice@epm.xsaic.com](mailto:mice@epm.xsaic.com).

If I can be of additional assistance on this and other matters, please write or call me at (202) 260-1639 at your convenience.

Sincerely,

Maria Gomez-Taylor, Ph.D.  
Analytical Method Staff  
Engineering and Analysis Division (4304)

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OCT 28 2003

OFFICE OF  
WATER

Les Orr  
Technical Representative  
Environmental Express  
490 Wando Park Blvd.  
Mt. Pleasant, SC 29464

Dear Mr. Orr:

This letter responds to your request for an evaluation of Environmental Express, Ltd.'s single use (disposable), carbon-coated polyethylene terephthalate (PET) BOD bottles. Generally, the ATP program evaluates only methods, not specific pieces of equipment. However, ambiguities in the approved methods made it unclear as to whether the single use, coated-PET bottles fell within the inherent flexibility of the methods. Therefore, EPA agreed to provide this evaluation.

Environmental Express provided data from seven laboratories comparing results obtained using single use, coated-PET bottles with those obtained using glass BOD bottles in a number of sample matrices. The data indicated that comparable performance was achieved with both bottle types. Therefore, EPA recommends that Environmental Express's single use, coated-PET BOD bottles be allowed for use with EPA-approved BOD methods (e.g., EPA Method 405.1) and associated methods for determination of dissolved oxygen (e.g., EPA Methods 360.1, 360.2) that are used in wastewater compliance monitoring. Each bottle should only be used for a single use, as directed by the manufacturer.

We appreciate your continued interest in environmental monitoring. If you have any questions regarding this letter or other issues, please contact me at 202-566061 at your convenience.

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style	Catalog number	size	reproducibility
ProWeigh	F93447MM	47mm	± 0.2mg
DoubleWeigh	F9344MM-X	47mm	± 0.2mg
ProWeigh for Volatiles	F93447VOL	47mm	± 0.2mg
ProWeigh	F93442MM	42mm	± 0.2mg
ProWeigh	F93455MM	55mm	± 0.2mg
ProWeigh	F93455VOL	55mm	± 0.2mg
ProWeigh	F93476MM	76mm	± 0.2mg
DoubleWeigh	F93476MM-X	76mm	± 0.2mg
ProWeigh for Volatiles	F93476VOL	76mm	± 0.2mg
ProWeigh	F93490MM	90mm	± 0.2mg
DoubleWeigh	F93490MM-X	90mm	± 0.2mg
ProWeigh for Volatiles	F93490VOL	90mm	± 0.2mg
ProWeigh	F934110MM	110mm	± 0.2mg
ProWeigh for Volatiles	F934110VOL	110mm	± 0.2mg
ProWeigh PES Filters 0.45µm	F96047PES	47mm	± 0.2mg

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